# Motion Pro Y Serie

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

## TABLE OF CONTENTS


- 3. Product and Standard Accessories
- 4. Getting Started
- 5. Camera Connector Legend
- 6. Status LED Definitions and Network Configuration
- 7. Daisy-Chaining
- 8-10. Dimensions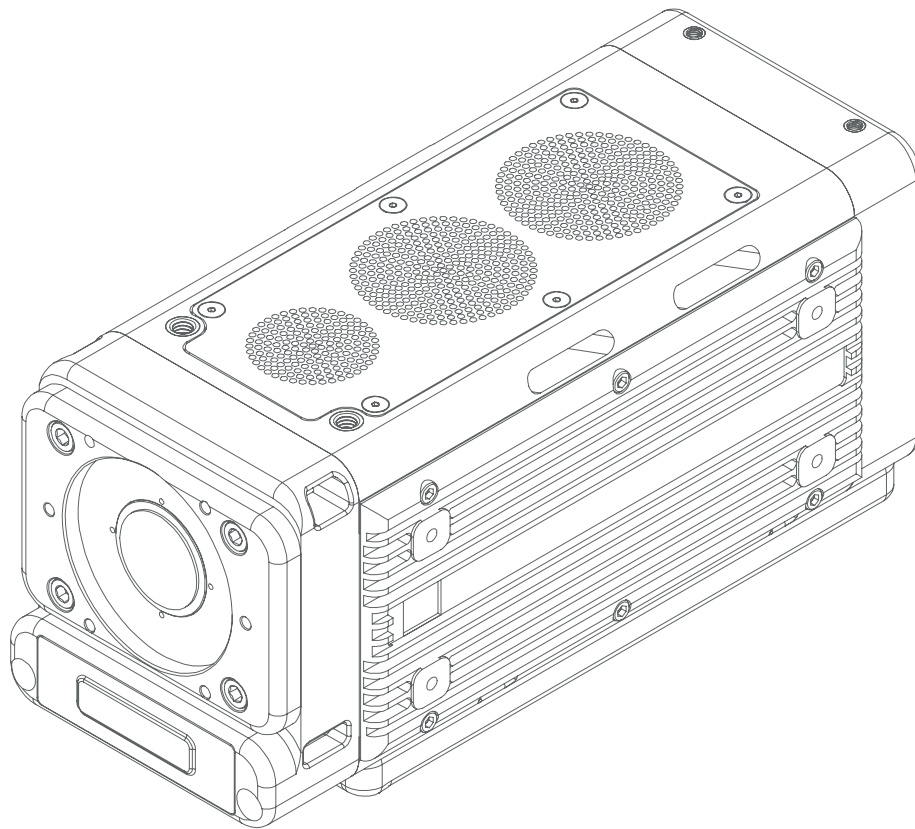


Y

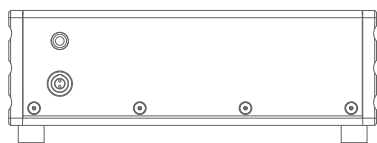Power Supply

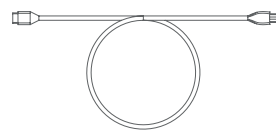
AC Power Cable

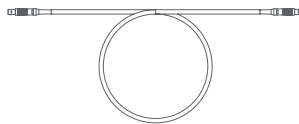
2-pin to 2-pin LEMO Cable

## Getting Started

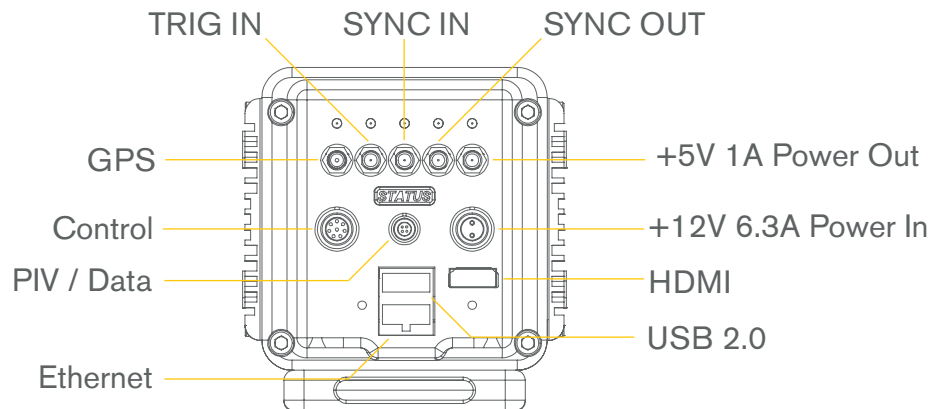CAUTION:
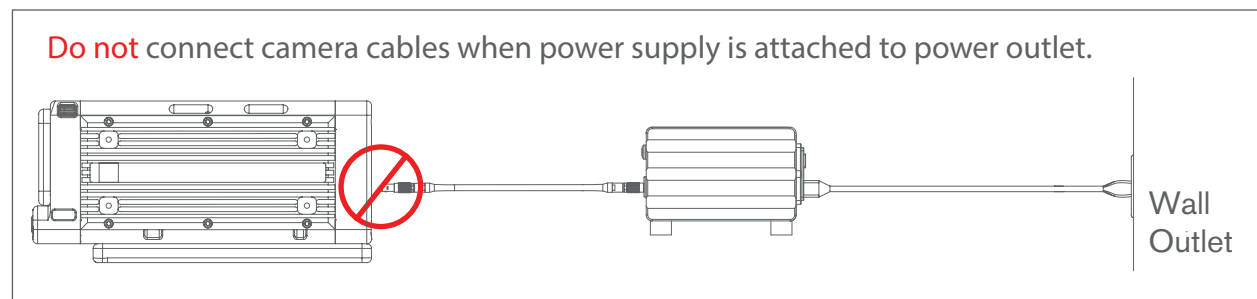
## Connecting the Y Camera to a PC or MAC computer:


**Caution:** Do not plug the Power Supply first. Be sure to follow the steps below.

**Note:** Descriptions for all connections can be found on the following page.

- Step 1** Make all SMA connections to the upper rear side of the Y Camera.
- Step 2** Connect Control, PIV, +12V 6.3A, HDMI, Ethernet or USB connections to the appropriate location on the rear of the Y Camera.
- Step 3** Using the 2-pin to 2-pin LEMO cable, line up the key of one LEMO connector and gently insert it into the 2-pin power connector on the back of the Y Camera until securely fastened.
- Step 4** Line up the key of the remaining LEMO connector to the 2-pin connector on the Power Supply and gently insert until securely fastened.
- Step 5** With the Power Supply switch in the 'OFF' position, connect the A/C Power Cable.
- Step 6** Once you have confirmed all connections are securely fastened, connect the AC Power cable plug in to a wall outlet.

# Camera Connector Legend

Definitions for the connectors on the back of the Y Camera:

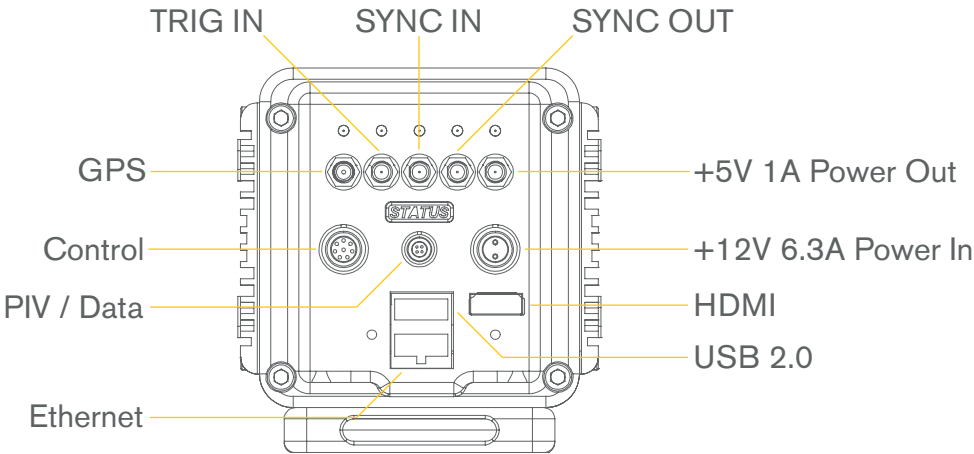**Trig in:** Connect an external Pickle Switch (5V TTL +/- signal or switch closure) to Trigger the camera to record images.

**Sync in:** Input signal, used for external synchronization.

**Sync Out:** Output signal, used for synchronization of multiple cameras in external sync mode or for synchronization with light source. Aligned with exposure time.

**GPS:** Connect an external GPS Antenna.

**Power Out:** +5V 1A SMA output, used for power of external WiFi module.

**Control:** Connector for IRIG (optional), Trig In, Sync In or Sync Out all differential.

**Power In:** Power input connector for the Y Series Power Supply.

**PIV / Data:** Connector for the use of Particle Image Velocimetry or external data peripheral device.

**HDMI:** HDMI output for an external display.

**Ethernet:** Ethernet connector to transfer data between the Y Camera and Computer and to control the camera.

**USB:** USB 2.0 connector to transfer data between the Y Camera and Computer and to control the camera.

## Status LED Definitions

The Y Camera has a status light on the rear of the camera. The following definitions indicate what status the cameras state is currently in:

When status LED color is mixed = Camera is booting up or processing

When status LED color is green = Camera is ready to record

When status LED color is red = Camera is recording images

## Network Configuration

**Note:** This information is for the use of the Ethernet connector on the Y Camera only.

Your Y Camera is delivered with an IP address schema as follows:

IP Address: 10.10.10.XXX

Subnet Mask: 255.255.255.0

**Note:** These properties can be changed but for the purposes of your initial configuration, you can leave them as is for now.

In order for your Y Camera to be detected by your computer, you need to confirm that the Network Card on your PC or MAC has an IP Address that falls within the range of the IP Address on the Y Camera. For example, if your Y Camera has an IP Address of 10.10.10.117, your computers network card needs to start with 10.10.10.XXX and have an ending numeric value from 1-255 but cannot be the same as the Y Camera IP Address (In this case the computers network card cannot be 10.10.10.117 as it will conflict with the IP address of the Y Camera).

**Note:** If using Ethernet, the network card on the computer that is connected to the Y Camera must support a Gigabit (1000BaseT) Ethernet connection.

## Daisy-Chaining

You can Daisy-chain multiple Y Cameras together in order to synchronize image recordings. This is useful if you want to film high speed footage from multiple angles simultaneously.


To Daisy-chain multiple Y Cameras together, make sure that the Y Cameras are powered down and follow these steps:

**Step 1:** Connect a SMA cable to the Sync Out connector on Y Camera 1.

**Step 2:** Connect the other end of the SMA cable to the Sync In connector on Y Camera 2.

**Note:** To Daisy-chain more than two cameras, simply repeat Step 1 and Step 2 above until you have connected all of the Y Cameras you wish to use.


Once you have completed all the steps above, you can follow the steps from page 4 to properly power up the Y Camera's.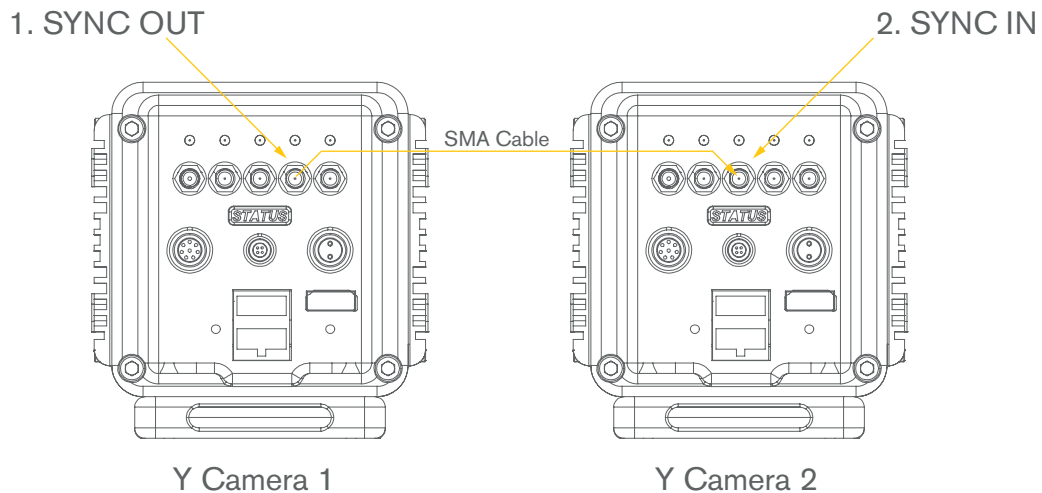

# Dimensions


## Front View

Unit of Measure =  $\frac{[mm]}{in}$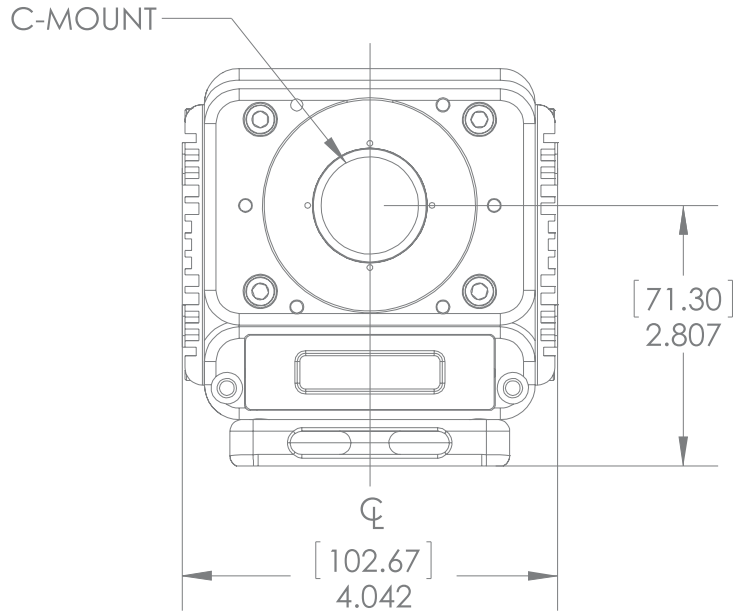

## Top View


Unit of Measure =  $\frac{[mm]}{in}$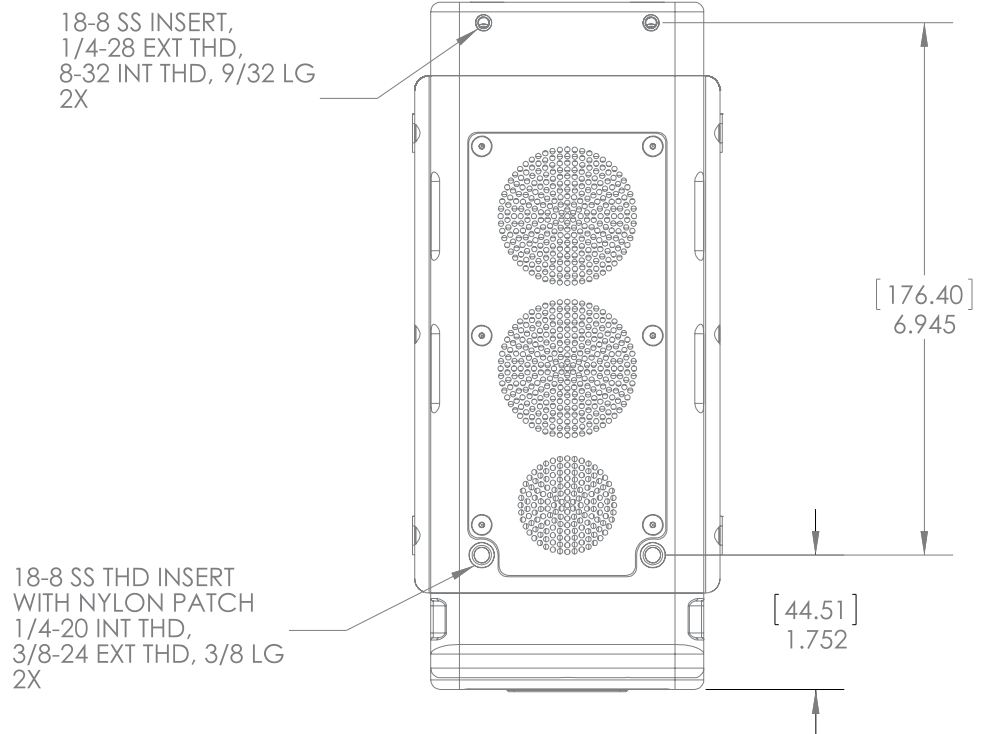

# Dimensions


Left Side View

Unit of Measure =  $\frac{[mm]}{in}$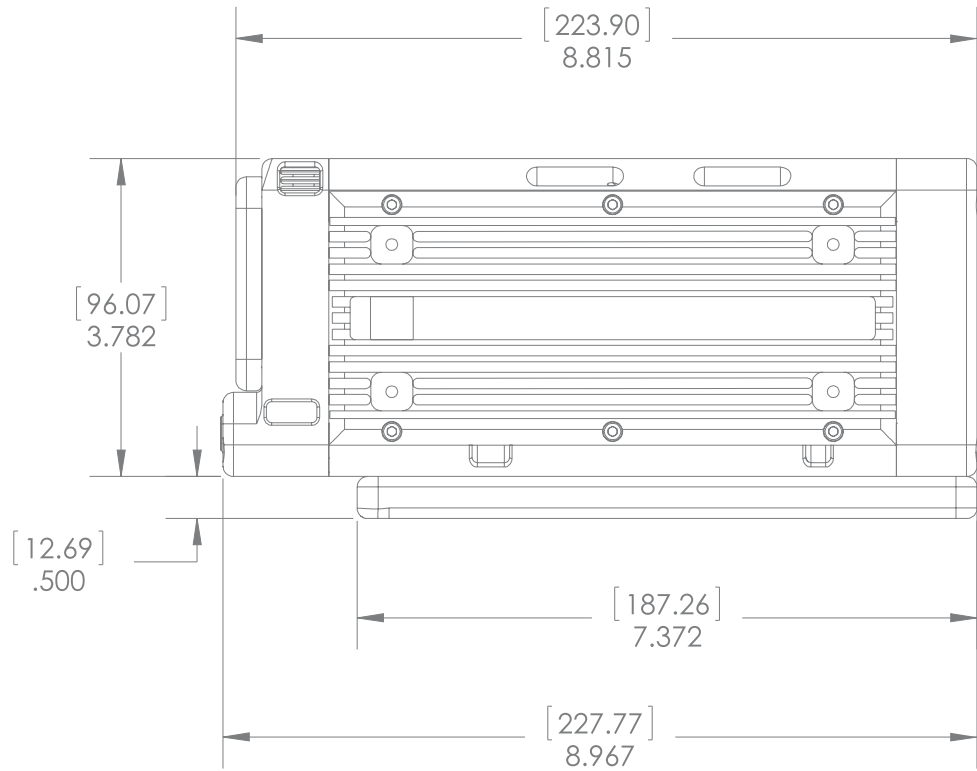

Right Side View


Unit of Measure =  $\frac{[mm]}{in}$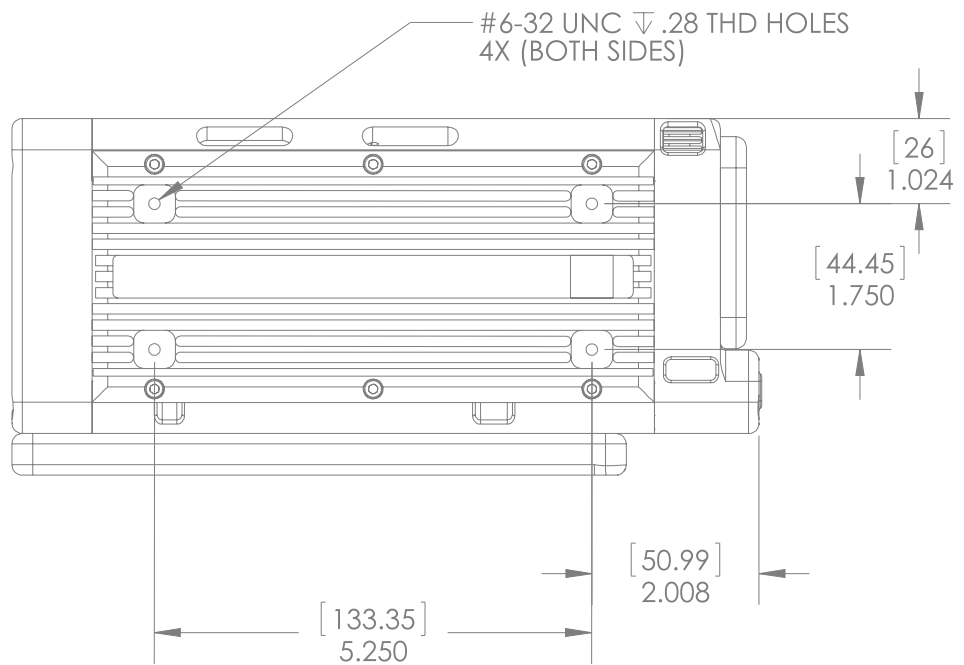

# Dimensions

## Rear View

Unit of Measure =  $\frac{[mm]}{in}$

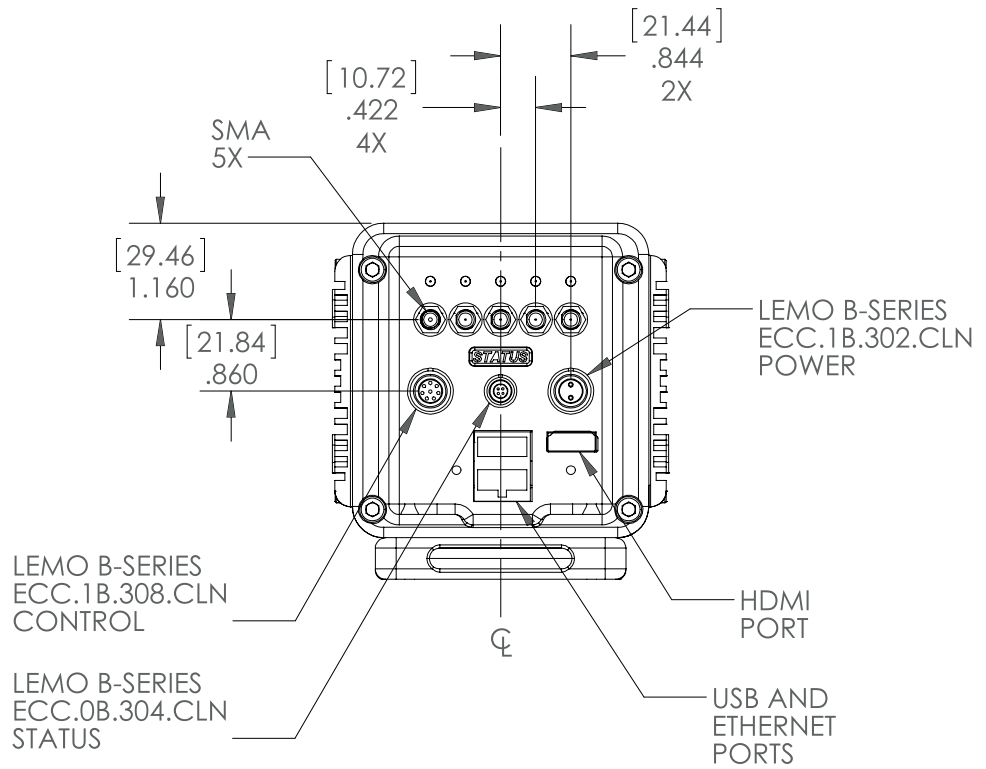## Bottom View

Unit of Measure =  $\frac{[mm]}{in}$